


PCR MYCOPLASMA DETECTION & ELIMINATION COMBO KIT

Mycoplasma detection is often difficult due to its lack of visible appearance, therefore it can be afflicting your valuable cell and affecting your results without your knowledge. Cells contaminated by mycoplasma species can have changes in proliferation, metabolism, gene synthesis and processing, and even adhesion properties. The solution for quick and easy mycoplasma detection is ABM's PCR Mycoplasma Detection Kit.

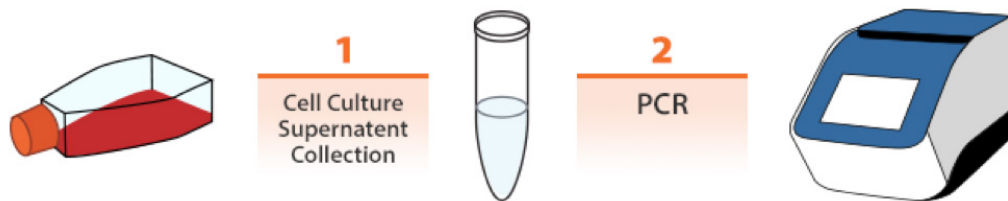

The **PCR Mycoplasma Detection Kit** allows for fast and reliable identification of mycoplasma contamination in cell cultures. Mycoplasma DNA in the cell culture supernatant is amplified via PCR and visualized using gel electrophoresis. In addition to the short detection process (less than 2 hours), the easy handling and high sensitivity makes this PCR Mycoplasma Detection Kit a convenient tool for routine examination of cell cultures and media.

PCR MYCOPLASMA DETECTION & ELIMINATION COMBO KIT	
Codice	Descrizione
G238	PCR Mycoplasma Detection Kit
G398	Mycoplasma Elimination Cocktail

Features

- Direct addition of cell culture supernatant to pcr reaction.
- No dna isolation/purification steps required.
- Ready-to-use primer mix.
- Reduces variability.
- Able to detect numerous mycoplasma species.
- High sensitivity.
- Included positive control to verify negative results.
- Rapid results in 2 hours.

Exclusive Distributor



Shipping Time
24h



Shipping Fees
€ 0



Shipping Fees for temperature controlled products
€ 0



Minimum Order
NO



La scelta logica

NAPOLI
Via Pisciarelli, 79
80078 - Pozzuoli (NA), Italia
Tel.: +39.081.610.7435
Fax: +39.081.610.7431

MILANO
Via Mauro Macchi, 8
20124 - Milano (MI), Italia
Tel.: +39.02.3823.4211
Fax +39.081.610.7431

offerte@microtech.eu
www.microtech.eu

PCR MYCOPLASMA DETECTION & ELIMINATION COMBO KIT

Number	Species and Strains
1	Mycoplasma verecundum
2	Mycoplasma verecundum strain ATCC 27862
3	Mycoplasma gallisepticum strain XZA
4	Mycoplasma gallisepticum strain JXB
5	Mycoplasma gallisepticum strain YMC
6	Mycoplasma gallisepticum strain ML
7	Mycoplasma gallisepticum strain WSA
8	Mycoplasma gallisepticum strain JXA
9	Mycoplasma gallisepticum strain WSB
10	Mycoplasma gallisepticum strain XZC
11	Mycoplasma gallisepticum strain F
12	Mycoplasma gallisepticum strain MG-HS
13	Mycoplasma gallisepticum strain YMA
14	Mycoplasma gallisepticum strain XA
15	Mycoplasma gypis
16	Mycoplasma gypis strain ATCC 51370
17	Mycoplasma indienne
18	Mycoplasma gallinaceum
19	Mycoplasma gallinaceum strain SA
20	Mycoplasma gallinaceum strain ATCC 33550
21	Mycoplasma gallinaceum strain 887
22	Mycoplasma gallinaceum strain 111R
23	Mycoplasma gallinaceum strain 33F
24	Mycoplasma arginini
25	Mycoplasma arginini strain EF-Hungary
26	Mycoplasma arginini strain ATCC 23243
27	Mycoplasma arginini clone V111-1
28	Mycoplasma sphenisci strain UCMJ
29	Mycoplasma sp. Mirounga ES2783_AG_GEN
30	Mycoplasma sp. Mirounga ES2774-NAS
31	Mycoplasma sp. Mirounga ES2806-NAS
32	Mycoplasma sp. Mirounga ES2774-NASSP4
33	Mycoplasma sp. Mirounga ES2714_AG_GEN
34	Mycoplasma sp. Mirounga ES2805-ORL
35	Mycoplasma pullorum
36	Mycoplasma pullorum strain ATCC 33553
37	Mycoplasma gallopavonis strain WR1
38	Mycoplasma gallopavonis strain ATCC 1197
39	Mycoplasma sp. CSL7489-oral
40	Mycoplasma sp. CSL7529-conjunctiva
41	Mycoplasma sp. CSL7491-lung
42	Mycoplasma sp. CSL7518-lung
43	Mycoplasma dispar strain ATCC 27140
44	Mycoplasma arthritis strain 91021
45	Mycoplasma arthritis strain ATCC 19611
46	Mycoplasma arthritis 158L3-1
47	Mycoplasma meleagridis
48	Mycoplasma meleagridis strain ATCC 27764
49	Mycoplasma iners strain PG30
50	Mycoplasma iners strain 89R-D-3
51	Mycoplasma iners strain 1D-E-3
52	Mycoplasma iners strain ATCC 15969

Number	Species and Strains
53	Mycoplasma fermentans M64
54	Mycoplasma fermentans JER
55	Mycoplasma fermentans strain ATCC 49892
56	Mycoplasma vulturni strain Gb-V33
57	Mycoplasma vulturni strain Gb-V33
58	Mycoplasma edwardii strain ATCC 23462
59	Mycoplasma edwardii strain 04-3436
60	Mycoplasma edwardii strain 04-3440
61	Mycoplasma flocculare
62	Mycoplasma flocculare strain Ms42
63	Mycoplasma conjunctivae HRC/581T
64	Mycoplasma conjunctivae strain Goat 655
65	Mycoplasma conjunctivae strain Goat 655
66	Mycoplasma conjunctivae strain HRC 581
67	Mycoplasma phocicerebrale ATCC 51405
68	Mycoplasma phocicerebrale strain 1049
69	Mycoplasma hyorhinis GDL-1
70	Mycoplasma hyorhinis clone V116-1
71	Mycoplasma hyorhinis strain GDL
72	Mycoplasma hyorhinis ATCC:29052
73	Mycoplasma hyorhinis isolate LuSIV
74	Mycoplasma hyorhinis strain ATCC 29052
75	Mycoplasma hyorhinis MCLD
76	Mycoplasma sp. Phocoena C-264-GEN
77	Mycoplasma sp. Phocoena C-269
78	Mycoplasma agalactiae
79	Mycoplasma agalactiae PG2
80	Mycoplasma agalactiae 5632
81	Mycoplasma neurolyticum strain ATCC 19988
82	Mycoplasma neurolyticum strain 10RN-F-3
83	Mycoplasma neurolyticum strain T15
84	Mycoplasma neurolyticum strain 36CFA2
85	Mycoplasma agassizii strain PS6
86	Mycoplasma agassizii strain Karina 09 Fr. Tortoise
87	Mycoplasma falconis ATCC 51372
88	Mycoplasma lagogenitalium strain ATCC 700289
89	Mycoplasma ovipneumoniae strain ATCC 29419
90	Mycoplasma columbinasale strain ATCC 33549
91	Mycoplasma felifaucium strain PU
92	Mycoplasma spermatophilum
93	Mycoplasma spermatophilum strain AH159
94	Mycoplasma spermatophilum strain 88(Jacobs)
95	Mycoplasma bovigenitalium strain MH 1113
96	Mycoplasma bovigenitalium strain PG-11
97	Mycoplasma bovigenitalium strain ATCC 14173
98	Mycoplasma bovigenitalium strain ATCC 19852
99	Mycoplasma equigenitalium strain ATCC 29869
100	Mycoplasma iguanae strain 2327
101	Mycoplasma testudineum strain ATCC 700618
102	Mycoplasma testudineum strain BH29
103	Mycoplasma moatsii strain MK405
104	Mycoplasma opalescens

Exclusive Distributor



Shipping
Time
24h



Shipping
Fees
€ 0



Shipping Fees for temperature
controlled products
€ 0



Minimum
Order
NO



La scelta logica

NAPOLI

Via Pisciarelli, 79
80078 - Pozzuoli (NA), Italia
Tel.: +39.081.610.7435
Fax: +39.081.610.7431

MILANO

Via Mauro Macchi, 8
20124 - Milano (MI), Italia
Tel.: +39.02.3823.4211
Fax +39.081.610.7431

offerte@microtech.eu
www.microtech.eu

PCR MYCOPLASMA DETECTION & ELIMINATION COMBO KIT

Number	Species and Strains
105	Mycoplasma zalophi strain CSL 4296
106	Mycoplasma zalophi strain CSL 5195
107	Mycoplasma zalophi strain CSL 4296
108	Mycoplasma gateae strain KDC
109	Mycoplasma gateae strain V-1
110	Mycoplasma gateae strain ATCC 23392
111	Mycoplasma canadense strain 466
112	Mycoplasma canadense strain 275C
113	Mycoplasma canadense strain ATCC 29418
114	Mycoplasma canadense strain 466
115	Mycoplasma corogypsi strain BV1-5
116	Mycoplasma corogypsi strain BV1-1
117	Mycoplasma equirhinis strain M432-72
118	Mycoplasma equirhinis strain ATCC 29420
119	Mycoplasma canis strain ATCC 19525
120	Mycoplasma canis strain C3b
121	Mycoplasma canis strain Skotti A
122	Mycoplasma hyopharyngis strain H36B-F
123	Mycoplasma hyopharyngis strain 538 Nasal
124	Mycoplasma auris
125	Mycoplasma cloacale
126	Mycoplasma mustelae
127	Mycoplasma mustelae strain MX9
128	Mycoplasma sp. 1220
129	Mycoplasma sp. 31848
130	Mycoplasma sp. 31948
131	Mycoplasma salivarium
132	Mycoplasma salivarium strain MP1166
133	Mycoplasma leonicaptivi strain ATCC 49890
134	Mycoplasma anatis
135	Mycoplasma anatis strain K6193C
136	Mycoplasma anatis strain K6193A
137	Mycoplasma felis strain ATCC 23391
138	Mycoplasma anseris
139	Mycoplasma hyopharyngis
140	Mycoplasma bovirhinis
141	Mycoplasma simbae ATCC 49888
142	Mycoplasma synoviae ATCC 25204
143	Mycoplasma synoviae 53
144	Mycoplasma orale
145	Mycoplasma orale strain MT-4
146	Mycoplasma hyosynoviae strain ATCC 25591
147	Mycoplasma bovoculi strain ATCC 29104
148	Mycoplasma maculosum strain ATCC 19327
149	Mycoplasma maculosum strain Skotti B
150	Mycoplasma elephantis strain ATCC 51980
151	Mycoplasma columbinum
152	Mycoplasma columbinum strain FG295
153	Mycoplasma caviae strain G122
154	Mycoplasma caviae strain ATCC 27108
155	Mycoplasma hominis ATCC 23114
156	Mycoplasma hominis strain ATCC 27545

Number	Species and Strains
157	Mycoplasma hyopneumoniae J
158	Mycoplasma hyopneumoniae 168
159	Mycoplasma hyopneumoniae strain ATCC 27714
160	Mycoplasma hyopneumoniae 7448
161	Mycoplasma hyopneumoniae 232
162	Mycoplasma zalophidermidis strain CSL 4779
163	Mycoplasma pulmonis strain Ash (PG34)
164	Mycoplasma pulmonis (strain UAB CTIP)
165	Mycoplasma alligatoris strain ATCC 700619
166	Mycoplasma crocodyli
167	Mycoplasma crocodyli MP145
168	Mycoplasma crocodyli strain ATCC 51981
169	Mycoplasma subdolium strain ATCC 29870
170	Mycoplasma phocicerebrale strain CSL 519552
171	Mycoplasma phocicerebrale strain ATCC 51405
172	Mycoplasma phocicerebrale strain Cheryl
173	Mycoplasma bovis
174	Mycoplasma bovis strain HEK-FDA
175	Mycoplasma bovis HB0801
176	Mycoplasma bovis PG45 clone MU clone A2
177	Mycoplasma bovis strain Madison
178	Mycoplasma bovis Hubei-1
179	Mycoplasma bovis strain 70-213 (Del Giudice)
180	Mycoplasma bovis strain ATCC 25025
181	Mycoplasma gallinarum strain ATCC 15319
182	Mycoplasma gallinarum strain B2 (Dinter)
183	Mycoplasma gallinarum ATCC 15319
184	Mycoplasma phocirhinis strain CSL7520-lung
185	Mycoplasma phocirhinis strain CSL 7475-4
186	Mycoplasma phocirhinis strain ATCC 49639
187	Mycoplasma collis strain ATCC 35278
188	Mycoplasma californicum strain ST6
189	Mycoplasma californicum strain ATCC 33461
190	Mycoplasma spumans
191	Mycoplasma buteonis strain Bb/T2g
192	Mycoplasma oxoniensis
193	Mycoplasma leopharyngis
194	Mycoplasma primatum strain HRC292
195	Mycoplasma primatum strain I30
196	Mycoplasma primatum strain ATCC 25948
197	Mycoplasma molare strain ATCC 27746
198	Mycoplasma lipofaciens strain ATCC 35015
199	Mycoplasma cricetuli
200	Mycoplasma glycyphilum
201	Mycoplasma citelli strain RG-2C
202	Mycoplasma citelli strain RG1D-L
203	Mycoplasma citelli strain RG1D-S
204	Mycoplasma citelli strain RG2C
205	Mycoplasma citelli strain RG1D-L
206	Mycoplasma columborale strain ATCC 29258
207	Mycoplasma columborale strain 18303
208	Mycoplasma sturni ATCC 51945

Exclusive Distributor



Shipping
Time
24h



Shipping
Fees
€ 0



Shipping Fees for temperature
controlled products
€ 0



Minimum
Order
NO



La scelta logica

NAPOLI

Via Pisciarelli, 79
80078 - Pozzuoli (NA), Italia
Tel.: +39.081.610.7435
Fax: +39.081.610.7431

MILANO

Via Mauro Macchi, 8
20124 - Milano (MI), Italia
Tel.: +39.02.3823.4211
Fax +39.081.610.7431

offerte@microtech.eu
www.microtech.eu